

MET/CAL 11.0.2 HOTFIX2 READ ME

May 30, 2024

This document includes instructions on applying HotFix2 to a MET/CAL Runtime and MET/CAL Editor version 11.0.2 deployment as well as a list of modifications included in this HotFix. This HotFix is cumulative and contains the changes made in the previous HotFix for MET/CAL 11.0.2.

Note: This HotFix only applies to MET/CAL version 11.0.2! Do not attempt to apply this HotFix to any other version of MET/CAL!

APPLYING HOTFIX2 TO MET/CAL 11.0.2

Preparing to apply the HotFix:

- 1) Download the MET/CAL 11.0.2 HotFix2 ZIP archive to a temporary folder.
- 2) Right-click the ZIP archive and select **Properties** from the popup menu. If there is a checkbox labeled **Unblock** on the lower right of the General tab, check this option and click OK. Otherwise, just click OK.
- 3) Extract the contents of the ZIP archive to a local folder.

To apply this HotFix to an existing MET/CAL workstation version 11.0.2:

- 1) Make sure MET/CAL Runtime and MET/CAL Editor are not running.
- 2) Copy the **mc.dll** file extracted from the archive to the following folders, replacing the existing files:
 - a. **C:\Program Files (x86)\Fluke\METCAL**
 - b. **C:\Program Files (x86)\Fluke\METCAL Editor\Extensions\METCAL.Editor.Project**

Repeat this process on all MET/CAL workstations when MET/CAL is installed.

MET/CAL 11.0.2 CUMULATIVE HOTFIX CHANGES

Key	Issue Type	Description
MC-13865	FIX	Resolved an issue with legacy calibrator FSCs when procedure is compiled that allow MOD1 to be taken from MEM.
MC-13845	FIX	Resolved an issue that prevented 5540A/5550A/5560A statements that included a coil from working properly.