

# MET/CONNECT 2.2.1 README

This document provides information for the MET/CONNECT 2.2.1 release.

## WHAT'S NEW IN VERSION 2.2.1?

- Added support for MET/CAL 11.2.0
- For a complete list, see the table below

## REQUIREMENTS AND CONFIGURATION DETAILS

The MET/CONNECT Server installation has the following requirements:

- MET/CONNECT version 2.2.1 can be deployed as a new system and can also update any prior MET/CONNECT version 1.x or 2.x system
- MET/CONNECT Server can only be installed on **64-bit Windows 10** or **Windows Server 2012 R2** or later operating systems
- MET/CONNECT Server requires an instance of **SQL Server 2016** or later. You must provide the installer with credentials to set up the database
- MET/CONNECT Server requires **Internet Information Services (IIS)**. The installer will attempt to enable and configure IIS automatically
- MET/CONNECT Server sets up a web site on the server for MET/CONNECT Service (database access for MET/CAL)
- MET/CONNECT Server sets up a web site on the server for MET/CONNECT API (for CMMS integration)
- MET/CONNECT Server sets up TCP port exception rules in the Windows Firewall for the MET/CONNECT Service (port 6382 or 44382) and MET/CONNECT API (port 6381)
- MET/CONNECT Server sets up a shared folder for and deploys MET/CAL shared files (\metcal)
- MET/CONNECT Server sets up a shared folder for MET/CONNECT files (\metconnect) including the MET/CAL Client installer
- In order to use SAML2 authentication, MET/CONNECT Service must be deployed as a secure web site using HTTPS

## MET/CONNECT 2.2.1 CHANGES

Key	Type	Description
METCON-3904	ADDED	Display a warning if the MET/CONNECT database is not at the expected version.

Key	Type	Description
MTM-13135	ADDED	Updated web.config file to ensure handler for *.svc is installed for new installations.
MC-14107	FIXED	Resolved an issue when using MET/CONNECT that would cause MET/CAL to enter a "not responding" state for a short period of time after aborting a procedure run before actual procedure execution begins.